SPECIFICATIONS Display: 3/4"-High 4-digit LCD

Timing Capacity: 99 min, 59 sec 1 second

0.01%

Press MIN and SEC buttons simultaneously to clear the

display to zero. The display may be cleared to zero

Accuracy: 30 seconds (audible) Alarm: Size/Weight: $2\frac{1}{2} \times 2\frac{1}{2} \times \frac{3}{4}$ ": 2 oz

CLEARING THE DISPLAY TO ZERO

whether or not the timer is running. TO SET TIMER

display.

Resolution:

 Press MIN, SEC buttons to set desired timing period. Press and hold MIN or SEC button to rapidly advance

TO START TIMER

- Once desired timing period is shown on display, press START/STOP button to start countdown (see Time-Out section).
- When timer reaches zero, the alarm sounds for 30 seconds then automatically shuts off unless turned off manually. Press any button to manually silence the alarm.

TIMEOUT

If a pause in the timing operation is required, press the START/STOP button to "freeze" the display at the current reading (time-out). To resume timing at the point it was stopped, press the START/STOP button again. Any number of time-outs may be performed.

BUILT-IN MEMORY FUNCTION

Pressing the START/STOP button to silence a sounding alarm also returns the display to the previously programmed time. Only the START/STOP button returns the display to the previously set time.

ALL OPERATIONAL DIFFICULTIES

If this timer does not function properly for any reason, please replace the battery with a new high quality battery (see Battery Replacement section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

A faint display, incorrect display, erratic display, or no display are all indicators that the battery must be replaced. On back of unit, slide off the battery cover. Remove the exhausted battery and replace with a new AAA alkaline battery. For proper placement, see drawing located inside battery compartment. Replace battery cover. Replacement battery: Control No. 1105.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

CONTROL COMPANY

12554 Old Galveston Rd. Suite B230 Webster, Texas 77598 USA Ph. 281 482-1714 • Fax 281 482-9448 E-mail sales@control3.com • www.traceable.com

Control Company is ISO 9001:2008 Quality-Certified by DNV and ISO/IEC 17025:2005 accredited as a Calibration Laboratory by A2LA.

TRACEABLE®
INSTANT-RECALL
MEMORY TIMER
INSTRUCTIONS

Traceable® is a registered trademark of Control Company. ©2016 (\$\frac{1}{4}\$) Control Company. 92-5122-00 Rev. 1 060816