***Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung***

FC4S - FC6S Flocculators

F105A0111, F105A0112

**General Information / Informazioni Generali / Informations Générales / Información General /
Allgemeine Hinweise**

Before using the unit, please read the following instruction manual carefully.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.

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This unit must be used for laboratory applications only.

The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions.

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This unit has been designed and manufactured in compliance with the following standards:

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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IEC/EN 61010-1

Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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During chemical-physical treatments of waters aimed to coagulate and precipitate pollutants, chemical coagulants are evaluated and their dosing optimized by jar tests. In relation to the variable composition of wastewaters the chemical dosing and the test frequency are to be adapted to local conditions. The jar test equipment manufactured by VELP is characterized by an electronic speed control with an independent speed settable for each place in order to optimize the result of settling so to reduce chemical consumption. FC4S and FC6S allow to pre-select 10 different speeds for each stirring point using the relative selectors located on the front panel (10-15-30-45-60-90-120-150-200-300). The steel structure epoxy painted is studied and tested in order to give the instrument an excellent resistance against the attack of chemical agents and corrosion in general. The stainless steel stirring rods are adjustable in height by a self-blocking chuck. The examined samples can be lighted by a back led lamp controlled by a switch located on front panel.

2. Assembly and installation

Upon receipt and after having removed the packaging, please check the integrity of the instrument. The box includes:

- Flocculator
- Plug for power supply
- Switching VELP 100-240V/12V without plug
- Instruction manual

2.1 Electrical connection

Before connecting the instrument to the power supply, make sure that the values on the rating plate correspond to those of the power supply. Connect the unit to the power supply using the switching.

Ensure that the socket and the relative cut-off device conform to current safety norms and easy to reach.

2.2 Start-up

Rotate the speed knob completely to the left. Place the beakers containing the sample and then, set the speed.

3. Operating controls

It's possible to select the desired stirrer speed by the knob "Stirrer rpm" on the frontal panel. It's possible to choose the desired values: 10-15-30-60-90-120-150-200-300 rpm by using the information written on the analogical scale around the knob.

At the end of the work cycle set the stirring speed knob on "0" rpm and turn off the light switch. If the instrument is not used for a long period it's better to disconnect the external power supply from the plug.

4. Maintenance

No routine or extraordinary maintenance is necessary apart from periodically cleaning of the unit as described in this manual. In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors. The instrument must be transported in its original packaging and any indications present on the original packaging must be followed (e.g. palletized).

4.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with an non-inflammable non-aggressive detergent.

5. Technical data

Power supply	100-240 V / 50-60 Hz
Power	18 W (FC4S) 23 W (FC6S)
Dimensions (WxHxD)	655x404x296 mm (FC4S) 935x404x296 mm (FC6S)
Weight (with plug)	12.5 Kg (FC4S) 18 Kg (FC6S)
Speed range	10-15-30-45-60-90-120-150-200-300 rpm
Construction material	Epoxy painted metal structure
Environmental ambient temperature	5 - 40 °C (41 - 104°F)
Storage temperature range	-10... + 60 °C (14 - 140°F)
Max humidity	80%
Pollution degree	2

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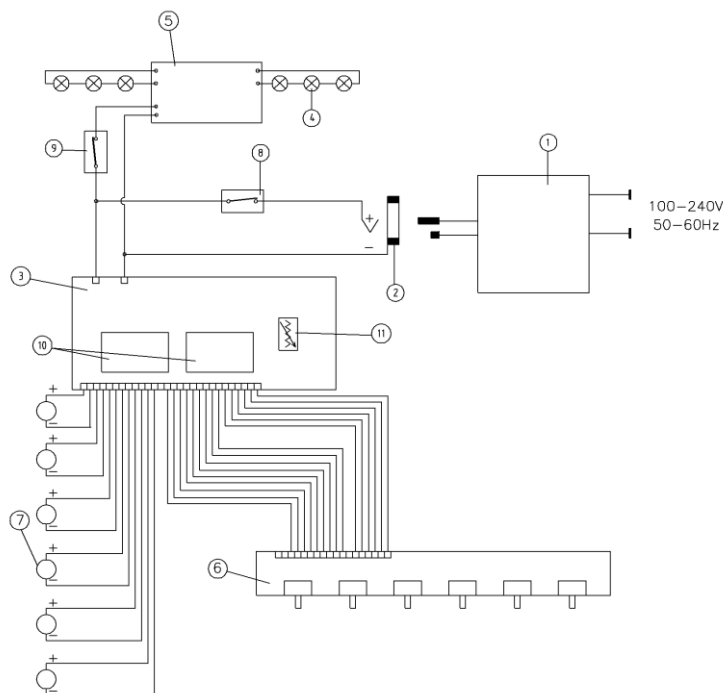
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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2011/65/EU (RoHS)

2012/19/UE (RAEE)

and satisfies the essential requirements of the following directives:

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- Machinery directive 2006/42/EC / Macchine 2006/42/CE / Machines 2006/42/CE / Máquinas 2006/42/CE / Maschinen 2006/42/EG

- Low voltage directive 2014/35/EU / Bassa tensione 2014/35/EU / Basse tension 2014/35/EU / Baja tensión 2014/35/EU / Niederspannung 2014/35/EU

- Electromagnetic compatibility directive 2014/30/EU / Compatibilità elettromagnetica 2014/30/EU / Compatibilité électromagnétique 2014/30/EU / Compatibilidad electromagnética 2014/30/EU / Elektromagnetische Verträglichkeit 2014/30/EU

- plus modifications / più modifiche / plus modifications / más sucesivas modificaciones / in der jeweils gültigen Fassung.

Thank you for having chosen VELP!

Since 1983 VELP has offered to professionals in the sector a range of sophisticated and reliable equipment.

VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Instruments are built according to the International norms IEC 1010-1 and to the rules of the CE mark.

Our product lines:

Food&Feed Line

DK and DKL Digesters
JP Recirculating water pump for fumes aspiration
SMS Scrubber
UDK Distillation units
SER Solvent extractors
FIWE Raw fiber extractors
Dietary fiber extractors
Oxitest Oxidation test reactor
NDA 701 Elemental analyzer
Consumables

Environment Line

ECO Thermoreactors
BOD. Determination systems
Refrigerated thermostats and incubators
Flocculators
Overhead mixer
Mineralization unit for trace heavy metals determination
Turbidimeter
Radiation detector

Stirring Line

Heating magnetic stirrers
Vertex digital thermoregulator
Ultraflat magnetic stirrer
Magnetic stirrers
Overhead stirrers
Heating plates
Vortex mixers
Homogenizer

Other Lab Solution

Open circulating bath
Recirculating water vacuum pump
Peristaltic pump

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