

I N S T R U C T I O N S



Replacement Hydrophobic Filters For Vacuum Aspirator Collection System

Not for use with HiFlow model, catalog no. F19917-0250

Catalog No. F19917-0155

These hydrophobic filters are to be used with the Scienceware® Vacuum Aspirator Collection System (Item # F19917-0150) in cases where collection bottle overflow has wet the filter and caused an interruption in vacuum. If you need instructions for the Vacuum Aspirator Collection System, please contact Bel-Art Customer Service at 1-800-423-5278 and ask for a copy of Instruction Sheet # 919917156.

NOTE: The hydrophobic filter protects the vacuum source from fluid contamination. Do not operate the pump without a hydrophobic filter in place as there will be no protection against fluids flowing into the vacuum pump and damaging the pump. Such damage is not covered by the warranty.

INSTRUCTIONS:

REMOVE OLD FILTER:

1. Grasp the filter and pull it upwards from the vacuum pump housing.
2. Unscrew the filter from the luer-lock contained within the tubing.
3. Discard the filter in accordance with all federal, state and local regulations.

INSTALL NEW FILTER:

4. Take one hydrophobic filter from the package and attach to the luer-lock by screwing the locking tabs on the filter into the luer-lock.
5. Press the other end of the hydrophobic filter into place in the small hole located on the vacuum pump housing above the On/Off switch.
6. You are now ready to resume normal operation of the Vacuum Aspirator Collection System.

If you need instructions for the Vacuum Aspirator Collection System, please contact Bel-Art Customer Service at 1-800-423-5278 and ask for a copy of Instruction Sheet # 919917156.

661 Route 23 South, Wayne, NJ 07470



TEL: 1-800-423-5278 • FAX: 973-694-7199 • www.belart.com

919917161 – 07/15

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products.

All such advice is given and accepted at the buyer's risk.

© 2015. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.

