




Item No.: H1000-MR

Flask Clamp Installation and Alignment


The MAGic Clamp Platform is capable of supporting a maximum of 5 x 500ml, 8 x 250ml, 12 x 125ml, or 20x50ml Erlenmeyer flasks. It is also capable of supporting various combinations of these four supported flask sizes.

To attach a flask clamp, simply place the clamp on the platform in the desired location. When attaching flask clamps for the maximum capacity of the platform, please follow the diagrams below for guidance:


5 x 500ml

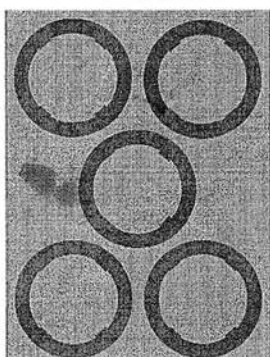


8 x 250ml

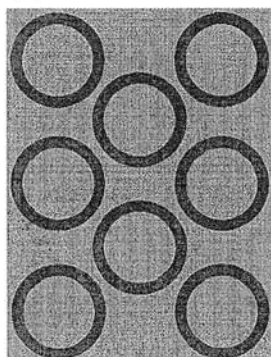

12 x 125ml

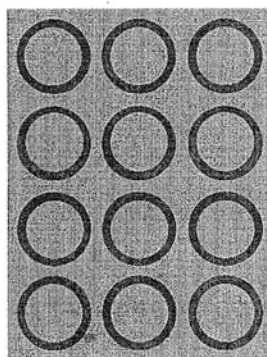

20 x 50ml

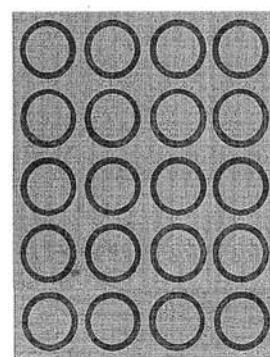

When attaching a mixture of different flask sizes, ALWAYS ensure a balanced load..

The MAGic Clamp platform is also capable of supporting various tube racks. The maximum capacity for each tube rack is listed below:

H1000-MR-1550	Combination rack for 15 & 50ml tubes	Max. 1
H1000-MR-0507	Tube rack for 72 x 5/7ml tubes (12-13mm)	Max. 2
H1000-MR-15ML	Tube Rack for 32 x 15ml tubes	Max. 2
H1000-MR-50ML	Tube Rack for 21 x 50ml tubes	Max. 2

