

Replacement of the Heating Element

1. Turn off BactiZapper unit and allow it to cool completely. Unplug the unit from the wall.
2. Remove the two screws holding the heating element to the base of the unit (Screws A):



3. Pull the unit off of the base gently, pulling up the electrical wires with it.
4. Unplug the head unit's two electrical wires from the base.
5. Plug the replacement element's two wires to those of the base and push the connected wires into the body of the BactiZapper until the replacement heating element is properly aligned with the body of the unit.
6. Replace the two screws removed in Step 2.