

LAB LABEL MAKER™ INSTRUCTIONS

Instructions: Loading tape– Set tape selector on side to 9.5mm; lift top black cover, insert tape cartridge colored side up, feed tape forward until tape is visible on white platform near the tip. *(May encounter some resistance, if difficulty feeding tape occurs, remove tape, curl from first 5 inches in the opposite direction to make tape straight).*

Printing tape– Slightly squeeze handle until it clicks to load tape and release handle; turn wheel to select character, completely and firmly squeeze handle; then select next character, to create a space or feed tape– slightly squeeze handle until it clicks to cut tape push firmly down on black lever at the very front lower tip of unit; peel off white backing.

Warranty, Service, or Recalibration

For warranty, service, or recalibration, contact:

CONTROL COMPANY

4455 Rex Road

Friendswood, Texas 77546 USA

Ph. 281-482-1714 Fax 281-482-9448

E-mail sales@control3.com

www.control3.com

Control Company is ISO 9001 Quality-Certified by DNV and ISO 17025 accredited as a Calibration Laboratory by A2LA.

Control Company Cat. No. 1600

Traceable® is a registered trademark of  Control Company

© 2009  Control Company 92-1600-00 Rev. 1 090809