

# Safety Pipette Fillers



A standard safety pipette filler eliminates the need for mouth pipetting

## Catalog No.

RED-001 Red Safety Pipette Filler  
BLK-001 Black Safety Pipette Filler  
BLU-001 Blue Safety Pipette Filler  
GRN-001 Green Safety Pipette Filler

## Features

100% Natural Rubber  
Standard Three Valve Balls  
1/4" Diameter Opening  
Up To 50ml Pipettes

- ❖ Also available in a silicone rubber version to provide greater chemical resilience and damage resistance
- ❖ Ask for catalog #SILRED for red or #SILBLK for black.

1. Squeeze the top valve (marked "A") and squeeze the bulb.
2. Let go of top valve before placing your pipette into the solution you intend to transfer.
3. Squeeze the "S" (middle valve) to pull the liquid into the pipette.
4. Squeeze the "S" (middle valve) when you want to empty the pipette while squeezing the bulb.
5. To force excess liquid from pipette, squeeze the "E" (side valve) while also squeezing the small bulb next to it.

