

## SPECIFICATIONS

<b>Timing Capacity:</b>	99 minute 59 second
<b>Alarm:</b>	60 Second beep duration
<b>Attachments:</b>	Stand/Magnet
<b>Functions:</b>	Count Down or Up
<b>Battery:</b>	1.5 Volt-LR44 Button Cell Auto Shut-off after 2 minute inactivity
<b>Memory:</b>	Retention During Shut Off (Push any button to turn on and return to previous setting)

## OPERATION

### 1. BATTERY INSTALLATION

Lay Timer face down and lift up stand. Slide open battery cover door. Remove the plastic battery

activation strip to activate battery. Digits on front of timer will appear

### 2. START/ STOP BUTTON

1. Press Start/Stop button to begin the countdown/count-up
2. Press Start/Stop again to stop countdown/count-up.

### 3. MIN/SEC BUTTON

Press MIN /SEC button to set minutes or seconds for countdown.

### 4. RESET TIMER

Press Min and Sec Buttons simultaneously to clear last setting memory and get back to zero for count up mode.

### 5. Fast Forward

Press and HOLD MIN or SEC button to fast forward numbers.


### ALL OPERATIONAL DIFFICULTIES

If this timer does not function properly for any reason, please replace battery with a new high-quality battery (see "Battery Replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing battery with a new fresh battery will solve most difficulties.

### BATTERY REPLACEMENT

An incorrect or weak display, no display, or operational difficulties indicate that the battery should be replaced. To open BATTERY COMPARTMENT,

slide BATTERY COMPARTMENT DOOR on back of unit in the direction of arrow. Remove exhausted battery and replace with any new 1.5 V button cell LR44 battery in appropriate positions to agree with polarity symbols (+ and -) in battery compartment. Replace battery cover.

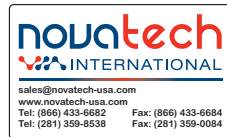

## WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

### TRACEABLE® PRODUCTS

12554 Old Galveston Rd. Suite B230

Webster, Texas 77598 USA

Ph. 281 482-1714 • Fax 281 482-9448

E-mail support@traceable.com • www.traceable.com

Traceable® Products is ISO 9001:2015 Quality-Certified by DNV and ISO/IEC 17025:2005 accredited as a Calibration Laboratory by A2LA.

Cat. No. 5128

Traceable® is a registered trademark of Cole-Parmer.

©2018 Traceable® Products. 92-5128-00 Rev. 0 011118

# TRACEABLE® BLUE-STEEL™ TIMER INSTRUCTIONS